

# Justifying a pathway to reduced replication-competent lentivirus (RCL) testing

*Adding the empirical to the theoretical.*

Vector manufacturing and analytics

14<sup>th</sup> May 2026

Dr. Dan Farley, Senior Director

Vector & Cell Innovation, OXB

Some graphics created in <https://BioRender.com>



# Legal disclaimer

This presentation does not constitute an offer to sell or a solicitation of offers to buy Ordinary Shares (the “Securities”). Although reasonable care has been taken to ensure that the facts stated in this presentation are accurate and that the opinions expressed are fair and reasonable, the contents of this presentation have not been formally verified by Oxford Biomedica plc (“OXB” or the “Company”) or any other person. Accordingly, no representation or warranty, expressed or implied, is made as to the fairness, accuracy, completeness or correctness of the information and opinions contained in this presentation, and no reliance should be placed on such information or opinions. Further, the information in this presentation is not complete and may be changed. Neither the Company nor any of its respective members, directors, officers or employees nor any other person accepts any liability whatsoever for any loss howsoever arising from any use of such information or opinions or otherwise arising in connection with this presentation.

This presentation may contain forward-looking statements that reflect the Company's current expectations regarding future events, its liquidity and results of operations and its future working capital requirements. Forward-looking statements involve risks and uncertainties. Actual events could differ materially from those projected herein and may depend on a number of factors.



# RCL testing impacts the full scope of LV product development

Why this talk has something for everyone



Review article

# A birdseye view of RCL testing

Will Development phase become a bottleneck?

0 1 2 3 4 (months to release)



Review article

Development phase

Qualification phase

Final RCL assay (routine testing)


Batch release



- **Amplification cell engineering?**
- Suitability/robustness?
- Sensitivity?
- Reg guidance



- Virus & cell banks
- Assay validation pack
- Inhibition studies
- Operator training
- SOPs & COPs

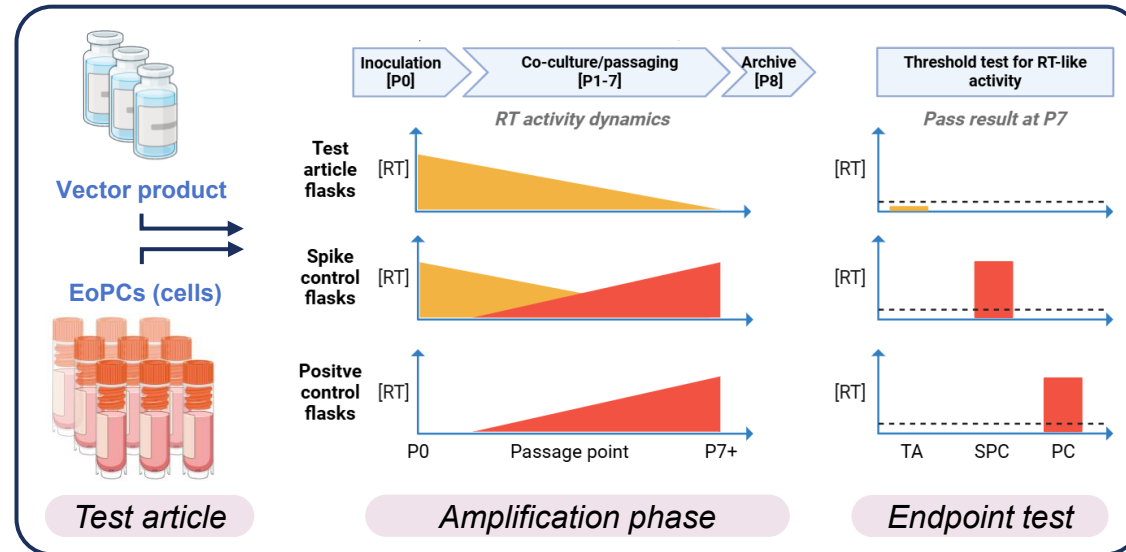


Test article

Amplification phase

Endpoint test

- Separate vector and EoPC tests
- Amplification phase =  $\uparrow$  RCL
- Positive & Spike controls
- Endpoint test on RT activity / p24 ELISA

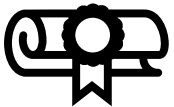

Pass result  
=  
Batch release

12+ months

6 months

2 months

➤ **A bespoke cell line for each retargeted LV envelope?** (RCL will most likely acquire envelope of the LV system employed)




# Safety features of contemporary LV systems

Why no RCL has been detected in their 25 year history



2KO/U1

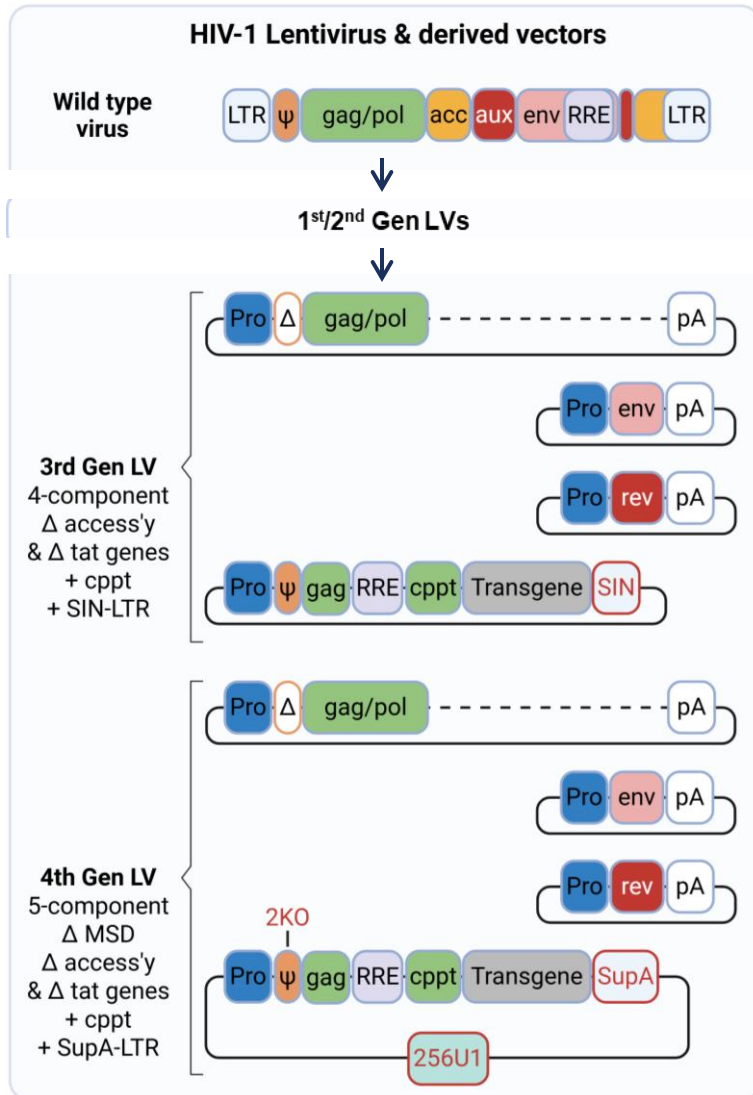

## Safer 3<sup>rd</sup> Gen RV/LV systems & improved in 4<sup>th</sup> Gen LV

- Components split into separate cassettes and plasmids
- Reduced homology (codon optimised GagPol – relevant later)
- Self-inactivating [SIN]-LTR & removal of tat
- Major splice donor mutation [2KO] and SupA-LTRs (OXB's SupA2KO-LVs)

➤ The more serial steps required to 'reverse' safety features, the less likely an RCL will form.




SupA-LTR





# A new model for assessing RCL formation risk

Possible vs Plausible vs Reasonable

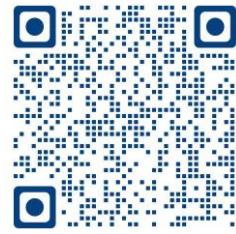

RCL model

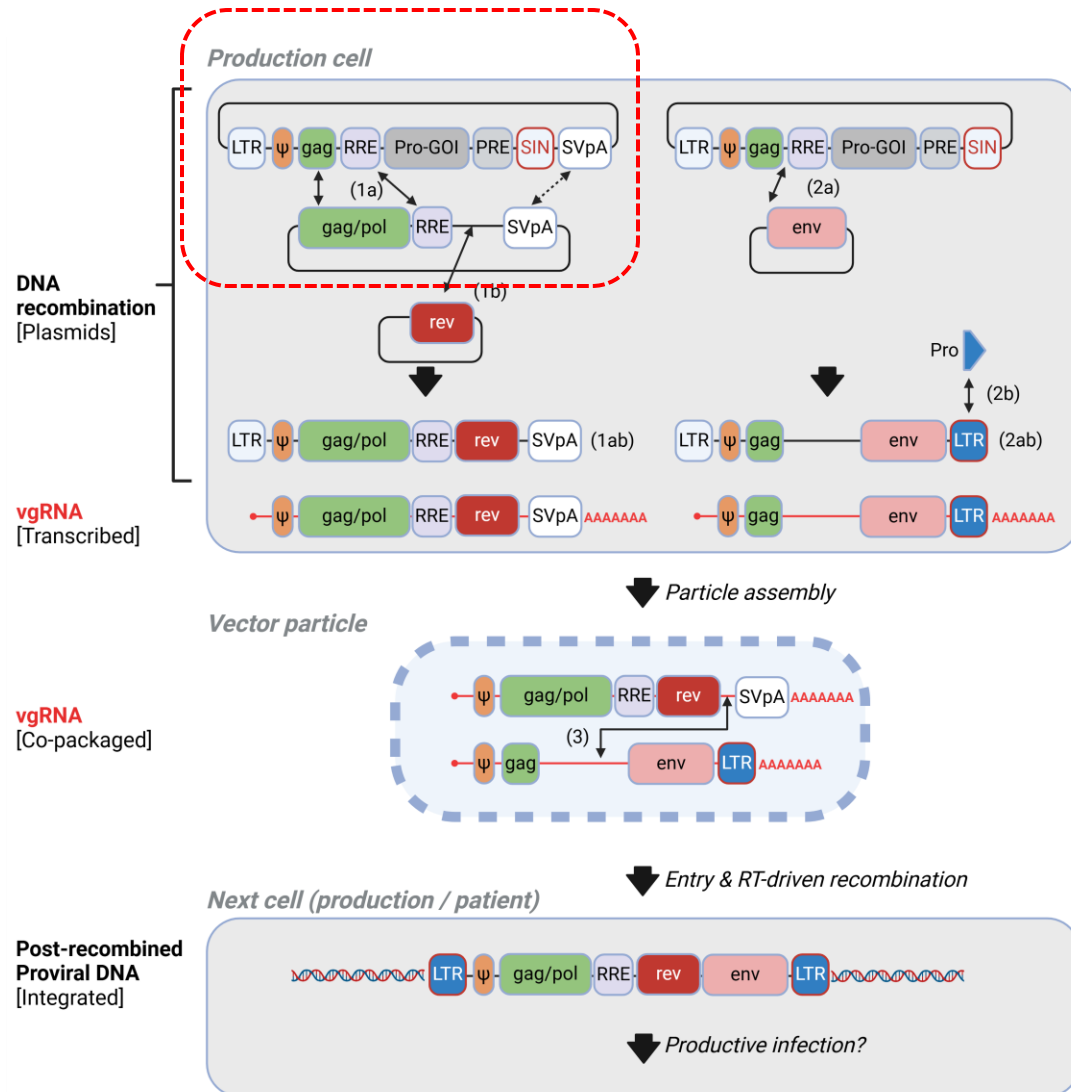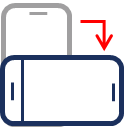


## The 'minimal' RCL model

- Probability of the arrival of a single RCL genome molecule in first cell
- Based on known cellular and HIV-1 biology mechanisms
- Using conservative assumptions and published data
- For 4x component transient:  $p[\text{RCL formation}] = 1$  in 10,000 bioreactors (200L)
- Probability reduces further with more components and more safety (SupA2KO-LV)
- **We are now adding empirical data from LV production cells, starting on Step 1a: genome x gagpol recombination....**

# Empirical detection of genome x gag/pol cassette recombination

A comparison of wild type HIV-1 vs codon-optimised gag/pol in HEK293T cells


## coGP has minimal homology to genome

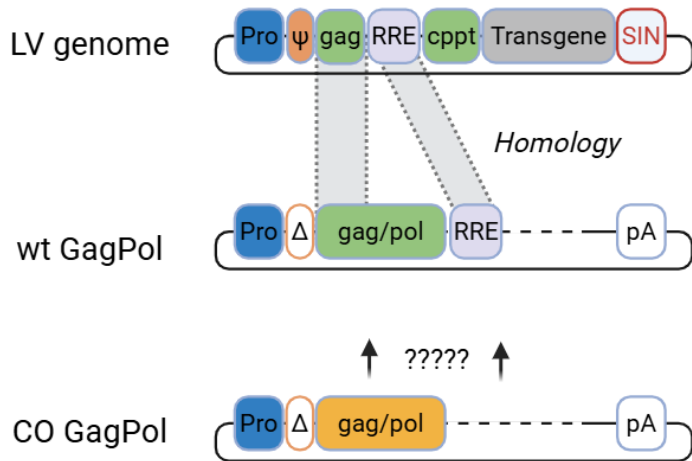

- wtGP shares Gag and RRE regions with genome
- coGP has minimal homology


- **Data indicates our predictions genome x wtGP recombinants per cell to be very close to reality (within 2-fold).**
- **Total pDNA is closer to our assumed transcriptionally active number (i.e. over-conservative)**

## LTR-Psi-Gag ddPCR product analysis

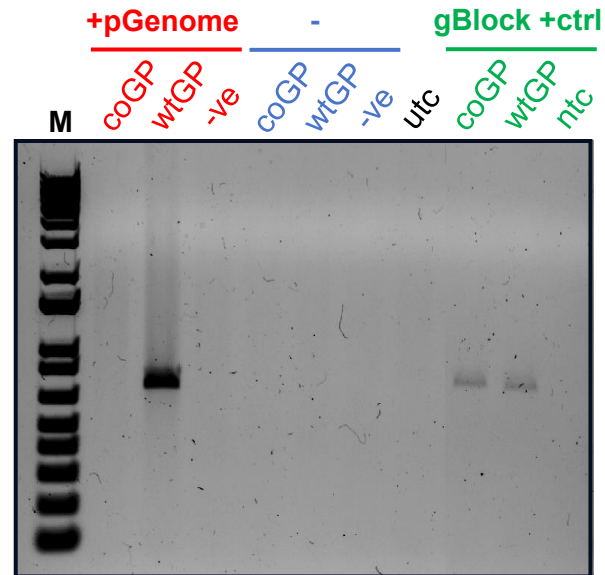

- Gel analysis of pooled ddPCR products
- Genome x wtGP recombination detected


## LTR-Psi-Gag ddPCR outcome

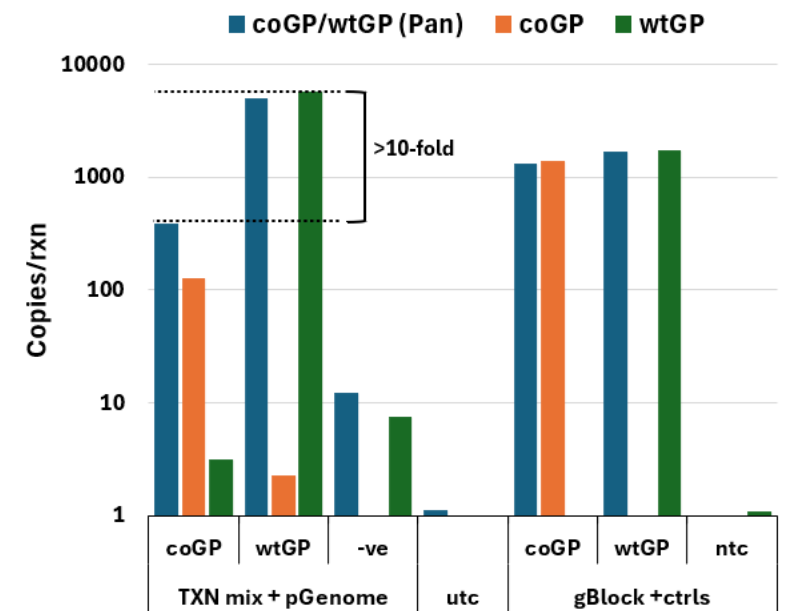

- ddPCR analysis show >10-fold lower signal for Genome x coGP
- Use of coGP strengthens case for reduced RCL testing

# Assessing risk from human endogenous RVs (hERVs)

Recombination risk from hERVs appears low

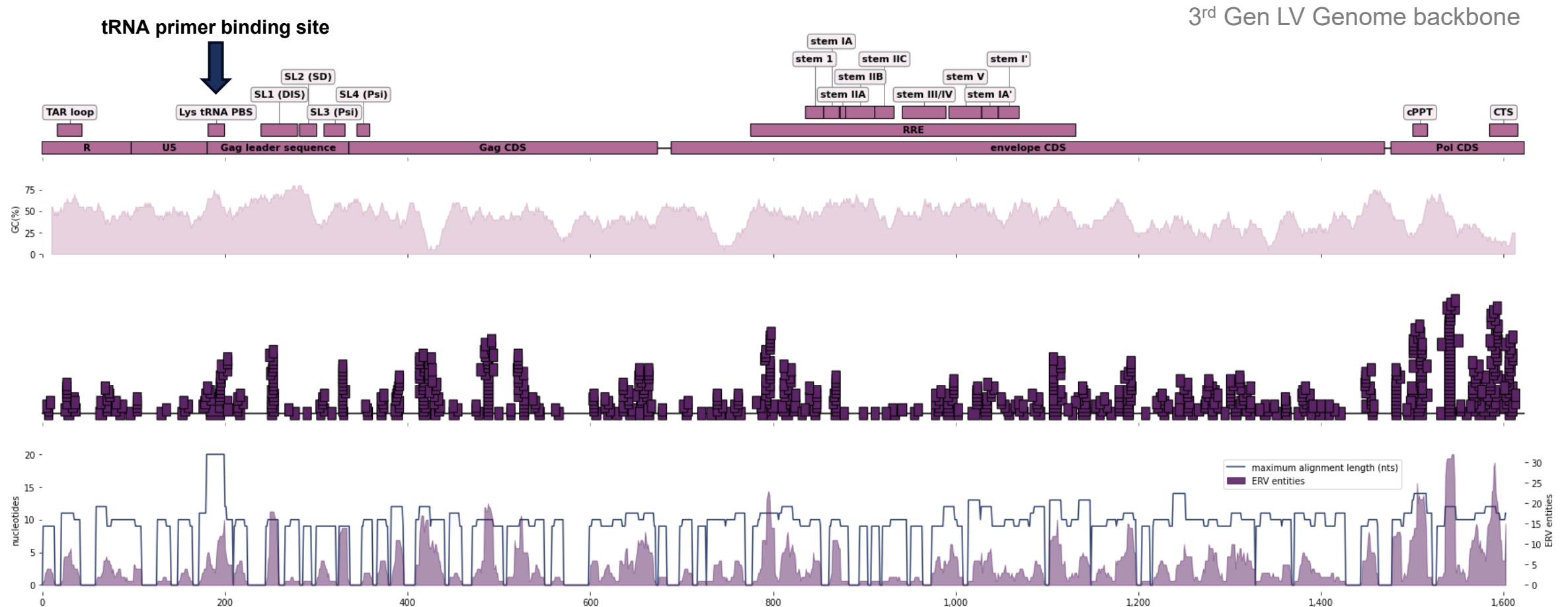


- Nucleotide alignment with **HERVd** database sequences (containing >0.5m endogenous HERV sequences) identified **HERV-K entities** as sharing greatest unbroken sequence homology with LV genome (~20 nts)
- Homologous alignments appear to cluster with **PBS**, **RRE** and **cPPT/CTS** regions

# Using the model to justify reduced RCL testing

A roadmap driven by data and probabilistic + logical arguments



Review article

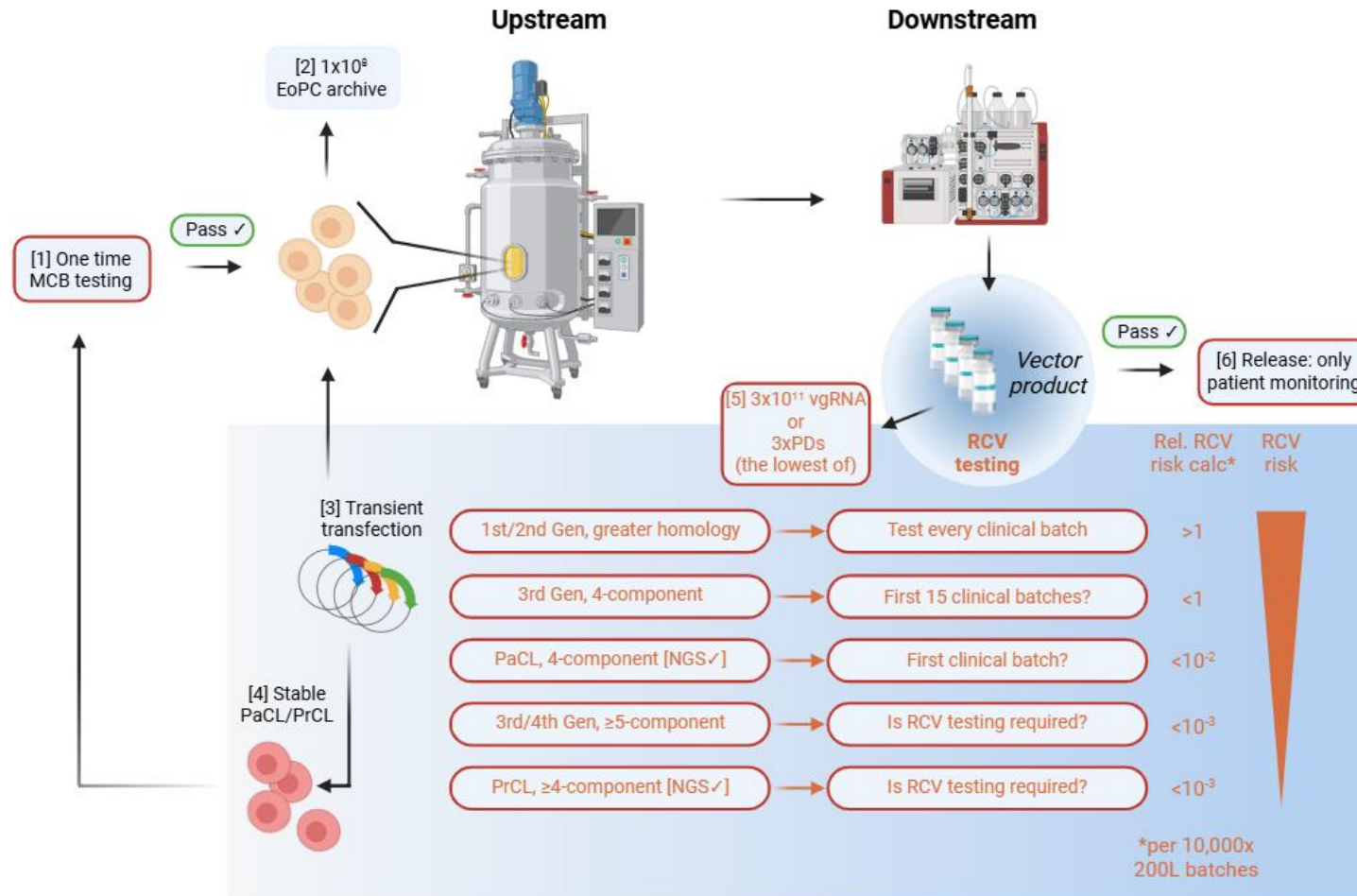

## Considerations for a roadmap to reduced / no testing

1. No more EoPC testing (archive cells only)
2. No testing of ex vivo modified cells (already close)
3. Fixed amount testing: 3-5x10<sup>11</sup> vgRNAs?
4. Fixed-and-finite testing? Only initial batches? 1-in-10 batches?
5. ≥5 components systems: only a supplementary assay?
6. PaCL and PrCLs: single, one-off testing?

# Launching the RADAR network

RCV Assay Development, Alignment & Regulation: consensus from the census



↑ Join RADAR ↑

1

## Connecting stakeholders from all parts of the industry

SMEs, Sponsors, CROs/CDMOs, Regulatory-advisors

2

## Surveys & feedback

Opportunity to feedback on the idea of the roadmap, RCL model, own experience and our surveys

3

## Sign-up for latest information

OXB feedback from surveys and progress

4

## Opportunity to input on navigating the regulatory pathway on RCL testing

If reduced testing is justified, how and who drives the change?

OXB

## OXB booth #1331

Come chat further, sign-up to RADAR, take the survey, get connected.



# Let's deliver life-changing therapies together

A global quality and innovation-led  
CDMO in cell and gene therapy

[OXB.com](https://www.oxb.com)



# Developing an RCL assay: no longer plug-and-play?


Beyond VSV-G; how bespoke, retargeted envelopes may bottleneck RCL testing




Review article

## Amplification cell and PC virus choices in Development

### LV/VSVG (off-the-shelf)

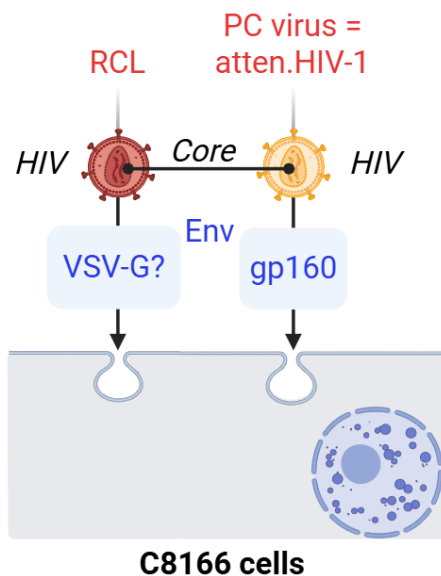


### Retargeted LVs

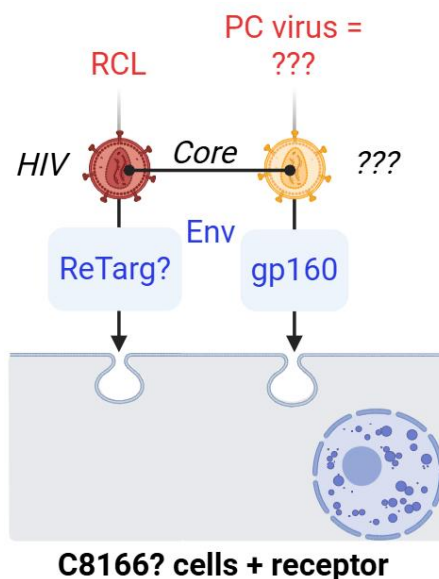

↑ **Essential properties** ↑

Desirable properties

## Bespoke RCL assay = Development bottleneck?

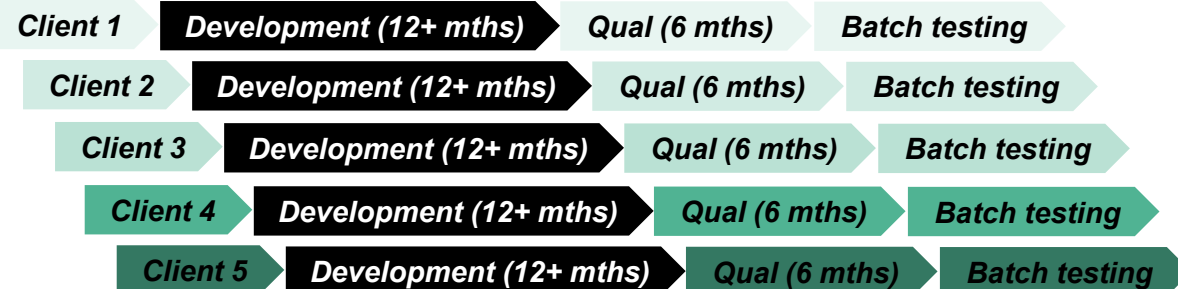

### Considerations for the CRO/CDMO, clients and the LV field regarding capacity


- Number project client batches?
- Number bespoke assays do you need to develop?
- The present solution is to amplify expensive RCL lab footprints
- Client expectation versus the reality of CRO/CDMO capacity.
- Competitive edge: non-viral platforms don't need any RCL testing!

➤ **Is this burden of testing justified?**



# What would we expect if an RCL was administered to a patient?

Reasonable inferences from what we actually know about HIV-1 and oncolytic RVs

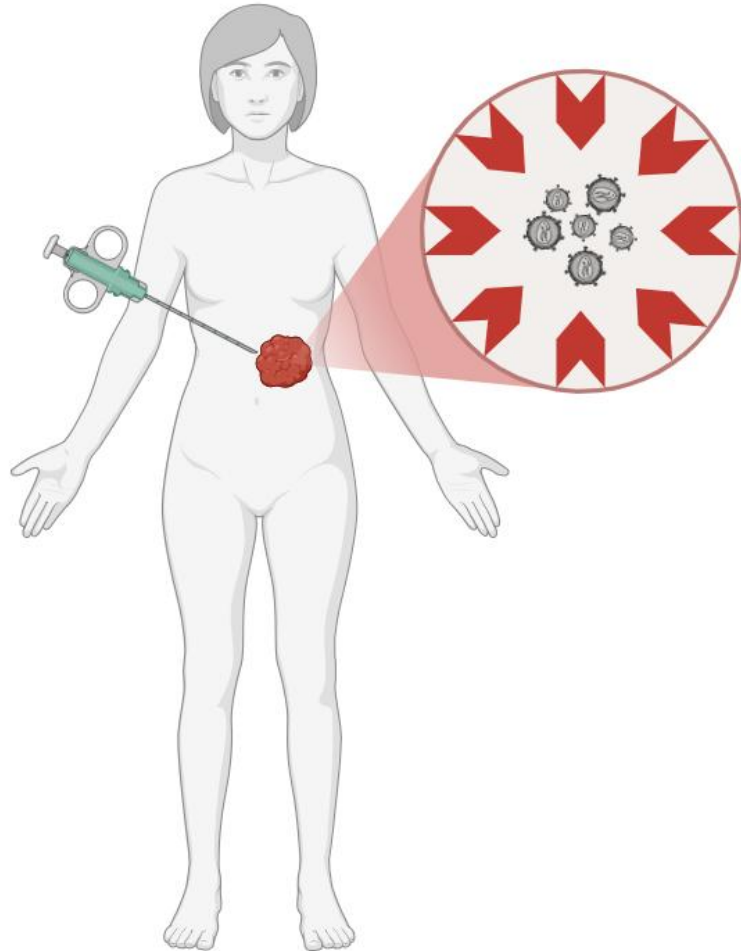

## Administration to patients

- **Highly unlikely** that a **minimal RCL** would be able to **replicate efficiently**
- Minimal RCL will **not encode accessory genes**: innate **restriction factors** will **block replication**
- No known strains of HIV-1 isolated from infected humans that **lack nef** lead to **pathogenic disease**
- **Vif and Vpr** expression are highly correlated with **lytic HIV-1 infection** of primary T-cells
- Minimal RCL will 'look' more like a simple RV like MLV: **MLV does not replicate in humans**
- MLV-based replicating **oncolytic RVs** (e.g. Toca 511) were injected into tumours (**no RCR testing was performed**, because they *are* replication-competent by definition!).
- Basis for **oncolytic RV safety** was in-patient 'containment' to tumour environment, relying on somatic tissue's **innate immune response** and slower cell division.